

BIO COOL HANDLING MANUAL

【Components Supplied】

- (1) EPS insulated shipper with air striation.
- (2) Plastic outer box or Carton outer box.
- (3) Labels includes: SHIPPING, UN3373, DRY ICE, OVER PACK.
(‘INFECTIOUS SUBSTANCE’ label is available. Please contact.)

【Packing Instruction】

Use for over pack of category A or third box of category B

- (1) Put the category A box or the secondary box of category B into the center of EPS inner box. Sometime you should fix these boxes to keep center.
- (2) Put dry ice, cool pack or heat storage pack into EPS inner box as necessary.
- (3) Close EPS top cover tightly and also carton.
- (4) Mark and label outer carton to applicable shipping regulations (IATA DGR).

【Label】

- (1) Fill in all the necessary items clearly on the ‘SHIPPING’ label, and put it on the corner of upper-left on carton.
- (2) Choose and put ‘INFECTIOUS’ or ‘UN3373’ label on the upper-right of outer carton.
In case of ‘UN3373’, put ‘BIOLOGICAL SUBSTANCE, CATEGORY B’ label close to the ‘UN3373’ label.
* Contact Specialists and get information about the classifications.
- (3) Put ‘DRY ICE’ label on the bottom-right corner if you need.
- (4) Put ‘OVER PACK’ label on the bottom-left corner.

【Safety precaution】

- (1) Do not use direct contact extreme cold like liquid nitrogen, high temperature water, chemicals and so on likely to give damages to EPS inner carton.
- (2) Prohibited the sterilization by ultraviolet and autoclave treatment.
- (3) Please use this BIO COOL box under specialists of transportation.
- (4) The classification of the specimen or the pathogenic agent should be advised by specialists.

Contact us: TREND SIGN CO., LTD.
6th floor, Nishi-Shinjuku TK Bldg., 8-19-2
Nishi-Shinjuku, Shinjuku-Ku, Tokyo 160-0023 JAPAN
TEL+813-5937-5721 FAX+813-5937-5722
info@bio-cool.net

Using the BIO COOL Box

(For Infectious Substance, category A and Biological Substance, category B)

